

Rapid Test Easy

High-performance qualitative assay



Cat.# M2241: Egg

Cat.# M2242: Casein

Cat.# M2243: Gluten

★Simple and Easy to use (Just 1-step)

★Specialized for swab and rinse water


★No need for special equipment (Swab kit is included.)

Performance	
Sample preparation time	Not needed
Assay time	10 min
Limit of detection	0.5 ppm
Test sticks	10 sticks/kit




Swab


Swab the specified area

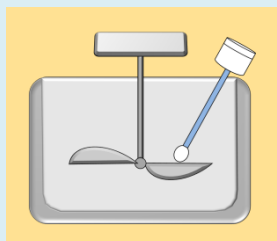

Shake vigorously

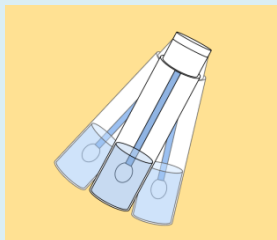


Add 200 μ L (4 drops) to the test stick

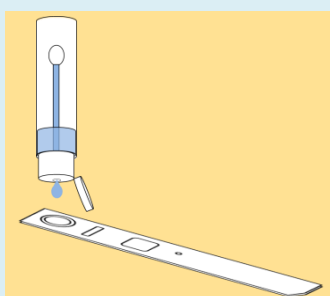


Rinse water

Collect rinse water

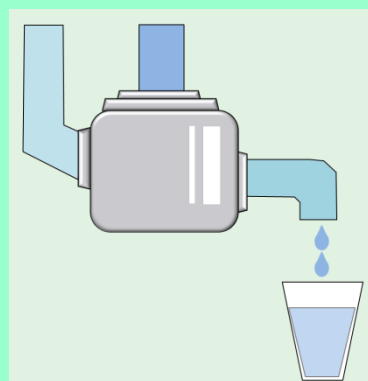

Simple, isn't it?

Add 200 μ L to the test stick

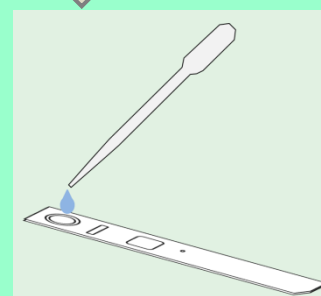

Just **10 min**
to check the result

Result

Positive

Negative



Component	Amount
Test stick	10 packs
Swab kit	10 bottles
Instruction	1 sheet