

Ultra Sensitive "PLUS" Mouse Insulin ELISA Kit



【Intended Use】

The Ultra Sensitive "PLUS" Mouse Insulin ELISA kit is for the quantitative determination of insulin in Mouse serum, plasma, and fluid. Please read the complete kit insert before performing this assay. The kit is for *RESEARCH USE ONLY*. It is not intended for use in clinical or diagnostic procedures or for internal or external use in humans or animals.

【Introduction】

Insulin is the primary hormone produced in the β cells of the Islets of Langerhans, and is known not only to regulate glucose metabolism, *i.e.* the uptake of blood glucose to the liver and peripheral tissues, but also play other important physiological roles.

Recent increases in the incidence of diabetes and obesity have stimulated intensive research on insulin levels and production. As a result, the accurate measurement of insulin in experimental animals is becoming increasingly important. The kit is a simple, precise, and sensitive ELISA sandwich assay for mouse insulin.

【Assay flow】

Preparation of standard and samples

↓
Dispense 45 μ L of Sample Diluent per well. Pipette 5 μ L of the sample (or working mouse insulin standard) per well.

First reaction

↓
Room Temperature, 2 hours.

Second reaction

↓
Room Temperature, 30minutes.

Enzyme reaction

↓
Room Temperature, 40minutes in dark.
Stop enzyme reaction.

Measurement of absorbance

【Precision】

The intra-assay precision (C.V.) : $\leq 10\%$

The inter-assay precision (C.V.) : $\leq 10\%$

The Lot to Lot precision (C.V.) : $\leq 10\%$

【Characteristics of Kit】

- **Sample size**
Only 5 μ L
- **Sample**
Serum, plasma, or fluid
- **Test**
96 wells (8 wells x 12 modules)
- **Reagents**
In liquid form (except standard)
- **Assay range**
0.025 - 1.6ng/mL
- **Assay time**
3hours 10min
- **Recovery**
When mouse insulin was spiked to mouse serum, the recovery of mouse insulin was 100% \pm 10%


【Specificity】

Substance	Reactivity
Mouse Insulin	100%
Rat Insulin	100%*
Hamster Insulin	Detected*
Rat C-peptide	Not Detected
Rat pancreatic polypeptide	Not Detected
Porcine glucagon	Not Detected
Human Insulin like growth factor-I	Not Detected
Human Insulin like growth factor-II	Not Detected

*Can vary from lot to lot.

Additional specificity data are available. In need, please contact to Morinaga Institute of Biological Science, Inc.

【Typical Standard】

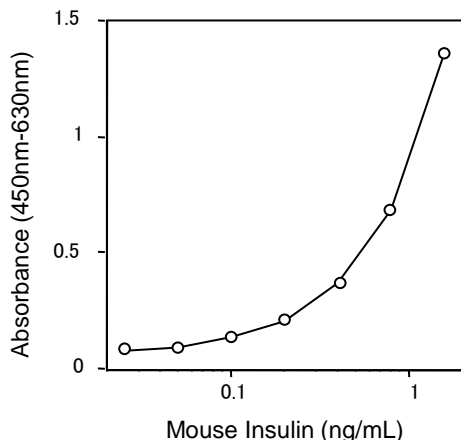

	Name of component	Content	Quantity
A	Antibody-coated microplate	96 wells	1 pack
B	Mouse Insulin Standard, Lyophilized	2.56ng/vial	1 vial
C	Standard Diluent	3 mL	1 bottle
D	Sample Diluent	13 mL	1 bottle
E	Anti-Insulin Enzyme Conjugate Solution	6 mL	1 bottle
F	Enzyme Substrate (TMB) Solution*	13 mL	1 bottle
G	Enzyme Reaction Stop Solution (1 N Sulfuric Acid)	13 mL	1 bottle
H	Wash Buffer Stock Solution (20X Concentrate)	50 mL	1 bottle
	Plastic microplate cover		1 piece

TMB solution* : Avoid exposure to light.

【Other required reagents and instruments】

- Micropipettes and disposable tips
- Volumetric flasks
- Distilled or deionized water
- Polypropylene microtubes
- Test tube racks
- Vortex mixer
- Aspirator for washing procedure
- Microplate reader
(capable of measuring A450 and A630 values)

【Contact us】

Morinaga Institute of Biological Science, Inc.
2-1-16 Sachiura, Kanazawa-ku
Yokohama 236-0003, Japan
e-mail: info@miobs.com