



V.1/Oct/2017

Quality Control for food allergen test

QC Material Egg
QC Material Milk
QC Material Wheat

***For Research or Laboratory Use Only.
Not for Use in Diagnostic Procedures.
Please read full descriptions in this instruction before use.***

Manufactured by:
Morinaga Institute of Biological Science, Inc. (MIoBS)
2-1-16 Sachiura, Kanazawa-ku, Yokohama 236-0003, Japan

E-MAIL : info@miobs.com
Website : <http://www.miobs-e.com>

I. INTENDED USE

The QC Material is the incurred food containing a certain amount of food allergen proteins. It can be used for skill evaluation, internal quality control, etc. of food allergen testing.

II. SPECIFICATION

1. This QC Material is a powder sample made of polysaccharide and sake residue.
2. The homogeneity is validated by the method referred to ISO13528.
3. The measurement values by the Food Allergen ELISA Kit II are shown on our web site (<http://www.miobs-e.com>).

III. LINE UP

QC Material Egg
QC Material Milk
QC Material Wheat

IV. COMPONENTS

Name of component	Content	Quantity
QC Material	Approx. 1.3g	5 packs
Instruction		1 copy

V. WARNINGS AND PRECAUTIONS

1. Use immediately after opening the package.
2. Treat the product with caution to avoid sample dispersion.

VI. HOW TO USE

1. Prior to use, bring the package to room temperature (20-30°C).
2. Weigh 1g each, and examine and evaluate with the method you execute routinely in the laboratory.

VII. STORAGE CONDITIONS AND EXPIRATION

1. Store at 2-8°C and avoid direct sunlight.
2. The expiration date (non-opened) is printed on the package.

VIII. WARRANTY

Morinaga Institute of Biological Science, Inc. makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. Buyer assumes all risk and liability resulting from the use of this product. There is no warranty of merchantability of the product, or of the fitness of the product for any purpose. Morinaga Institute of Biological Science, Inc. agrees to replace any defective product, but expressly disclaims liability for damages, including special or consequential damage, or expenses arising directly or indirectly from the use of this product.