

Rapid Test Easy for food allergens

High-performance qualitative assay



Cat.# M2241: Egg

Cat.# M2242: Casein

Cat.# M2243: Gluten

Cat.# M2244: Buckwheat

Cat.# M2245: Peanut


Cat.# M2246: Soya

Cat.# M2247: Crustacean

- ✓ **Simple and Easy to use** (Just 1-step)
- ✓ **Specialized for swab and rinse water**
- ✓ **No need for special equipment** (Swab kit is included.)


Performance

Sample preparation time	Not needed
Assay time	10 min
Limit of detection	0.5 ppm (μg protein /mL)
Test sticks	10 sticks/kit

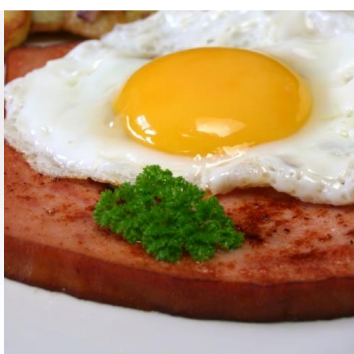


Swab


Swab the specified area

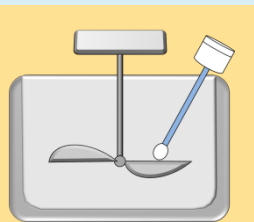

Shake vigorously

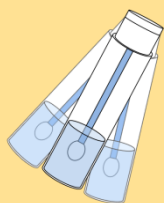


Add 200 μ L (4 drops) to the test stick

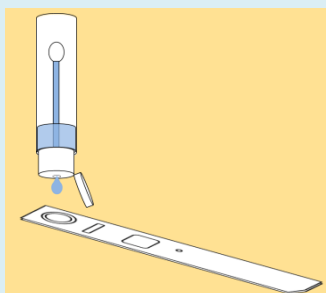


Rinse water

Collect rinse water

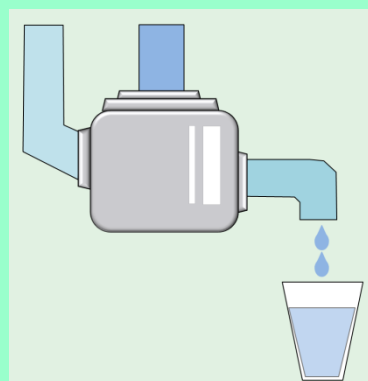

Simple, isn't it?

Add 200 μ L to the test stick

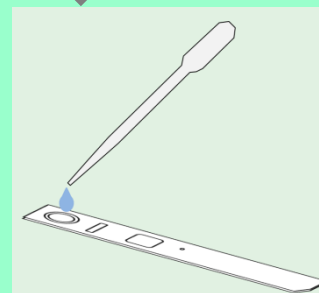

Just **10 min**
to check the result

Result

Positive

Negative



Component	Amount
Test stick	10 packs
Swab kit	10 bottles
Instruction	1 sheet