

Morinaga Food Allergen Lateral Flow  
Rapid Test Pro II for Shellfish

Rev. Sep. 2022

Lot number from #PL001

Group	Ingredient	Test Result	Group	Ingredient	Test Result	Group	Ingredient	Test Result	
<b>Cereals</b>	Rice powder	-	<b>Seeds</b>	Peanut	-	<b>Fruits</b>	Orange	-	
	Polished rice	-		Roasted almond	-		Kiwi	-	
	Sprouted brown rice	-		Roasted cashew nut	-		Apple	-	
	Glutinous rice	-		Roasted macadamia nut	-		Peach	-	
	Buckwheat	-		Roasted pistachio nut	-		Banana	-	
	Japanese barnyard millet	-		Roasted walnut	-		Coconut milk	-	
	Foxtail millet	-		Roasted pecan nut	-	Coffee bean	+		
	Common millet	-		White sesame	-	Cocoa powder	-		
	Wheat	-		Black sesame	-	Green tea	+		
	Rye	-		Salmon	-	Roasted laver	-		
	Barley	-	Mackerel	-	Hijiki (boiled)	-			
	Malt	-	Squid	+	Brown seaweed (raw)	-			
	Oat	-	Octopus	+	Kelp 200-fold extract	-			
	Adlay	- *	Black tiger prawn	+	Cumin	-			
	Corn	-	Snow crab	+	Poppy seeds	-			
	Corn flour	-	Scallop adductor	+	Black pepper	-			
	Amaranth	-	Pollack roe	-	White pepper	-			
Quinoa	-	Salmon roe	-	Red pepper	-				
<b>Tubers· Starch</b>	Sago palm powder	-	<b>Meats</b>	Beef	-	<b>Spices</b>	Ginger	-	
	Yam	-		Chicken	-		Cinnamon	-	
	Potato starch	-		Pork	-		Coriander	-	
<b>Beans</b>	Soybean	-	<b>Eggs</b>	Chicken egg	-		<b>Polysaccharide Thickener</b>	Turmeric	-
	Red bean (azuki)	-		Processed chicken egg	-			Low methoxyl pectin 50-fold extract	-
	Common bean	-		Dried chicken egg	-		Guar gum 1000-fold extract	-	
	Concord pole	-	<b>Dairy products</b>	Milk	-		<b>Others</b>	Gelatin	-
	Broad bean	-		Skim milk	-				
<b>Green stuff</b>	Tomato	-	<b>Fungus</b>	Mushroom	-				
	Onion	-							
	Spinach	-							

Test Result	
-	Negative
+	Positive

\*The Test Solution may be difficult to flow on a test stick.

These samples were tested by following the instruction of Rapid Test Pro II (using Recommended Extraction Method), except Kelp 200-fold extract, Polysaccharide and Thickener samples.

These results can be changed by the difference of kit's lot number or the sample's species or production area.